Readout

HORIBA Technical Reports

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January 1992 ■ No.4

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Why is Particle-Size Analysis so Difficult ?

There are many particle-size analyzers available on the commercial market and these various analyzers are based on a number of diverse principles. The situation tends to confuse potential users of particle-size analyzers : Why are there so many different kinds of equipment and what are the differences between them? It is also problematic that the various types of analyzers show discrepancies in the results of measurements.

One reason there are so many different types of equipment on the market is that the range of particle sizes that must be measured in today's particle-size analysis is simply too extensive to be handled by a single type of device. In addition to this, there are two reasons for discrepancy in results from device to device : (1) differences in the principles of measurement, particularly in the method of treating the optical properties of particles, and (2) differences in how the sample is prepared for measurement.

Of these three problems, one appears unavoidable : discrepancies resulting from different principles of measurement. By making efforts toward greater standardization, however, we should be able to overcome the other two problems, i.e., the various kinds of equipment on the market and the differences in the method of sample preparation.

Standardization is a goal we should be aiming at right now, and also, to ensure that we are responding adequately to the needs of potential users of our equipment, we should make sure that even if a great variety of analyzer equipment is to remain on the market, the principles and purposes of each type of device should be explained clearly.