

DSS-SPSE020

Two Color Silicon and Lead Selenide Solid State Detector

ELEMENTAL ANALYSIS
FLUORESCENCE
GRATINGS & OEM SPECTROMETERS
OPTICAL COMPONENTS
FORENSICS
PARTICLE CHARACTERIZATION
RAMAN
SPECTROSCOPIC ELLIPSOmetry
SPR IMAGING


An excellent two-color detector composed of silicon and lead selenide photodiodes, available in two different cooling levels, provides broad spectral response from the UV to near-IR.

These dual color single point detectors detect signals across a broad wavelength range, with a dual layered photodiode design using silicon on top of indium gallium arsenide. The silicon photodiode is transparent to wavelengths above 1000 nm allowing the InGaAs to detect those longer NIR wavelengths. With high sensitivity (D^*) and two options for ambient and thermoelectric cooling, responsivity extends from 200 nm to 2600 nm. This is one of a number of dual color single point solid state detectors available from HORIBA Scientific. Contact us for further information.

Used in conjunction with optically optimized housings, these detectors integrate seamlessly with HORIBA's extensive selection of monochromators. In addition, the SpectrAcq2 acquisition module allows for software integration with LabSpec, SynerJY, or LabVIEW. With all of the additional Optical Building Blocks available from HORIBA, a user can easily go from individual components to a complete spectroscopy solution.

Features and Benefits


- Wide spectral responsivity from 200 nm to 4500 nm
- High sensitivity ($D^* \sim 10^{12}$)
- Compact ambient and TE detector housing

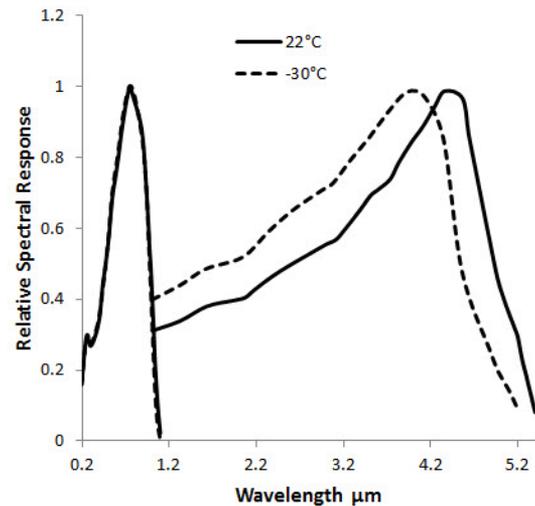

Accessories

Various accessories are available for powering the detectors, optically coupling detectors to HORIBA monochromators, and data acquisition.


- Power supply for TE cooled detector, DSS-15V-TEP
- Power supply for ambient, DSS-15VP
- Mirror-based housing, 1427C
- BNC cable, J30646
- SpectrAcq2 data acquisition module
- SMA fiber adapter, DSS-SMA
- Dual 1427C housing adapter, J23078370
- Dual detector housing, J23079050
- BNC switchbox for dual detectors, SWB-AB

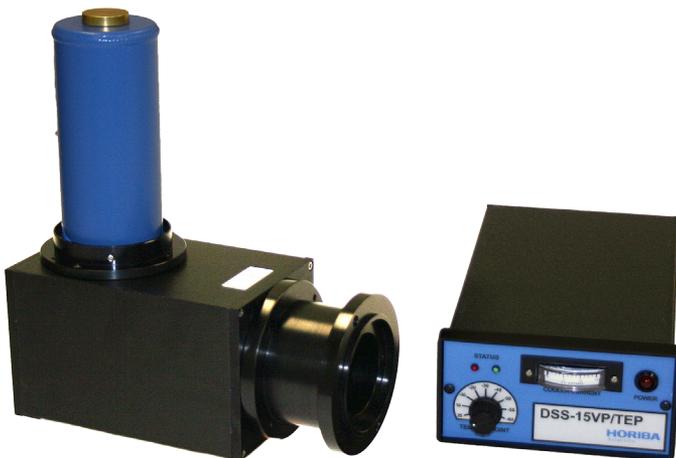


Specifications

Part number	DSS-SPSE020A		DSS-SPSE020T	
Detector type	2.5 mm diameter silicon photodiode / 2 mm x 2 mm lead selenide photoconductor			
Operating temperature (°C)	22°C ambient		-30°C TE cooled	
Operating wavelength (µm)	0.2 – 1.0 µm	1.0 – 4.5 µm	0.2 – 1.0 µm	1.2 – 4.7 µm
Responsivity (V/W @ peak)	0.6 x 10 ⁸ / 10 ⁷	1 x 10 ⁵ / 10 ⁴	0.55 x 10 ⁹ / 10 ⁸	2 x 10 ⁷ / 10 ⁶
Noise (V/Hz^{1/2})	1.2 x 10 ⁻⁶ / 10 ⁻⁷	10 x 10 ⁻⁶ / 10 ⁻⁷	3.6 x 10 ⁻⁶ / 10 ⁻⁷	10 x 10 ⁻⁶ / 10 ⁻⁷
NEP pk, (W/Hz^{1/2})	< 2.0 x 10 ⁻¹⁴	1.0 x 10 ⁻¹⁰	6 x 10 ⁻¹⁵	1 x 10 ⁻¹²
Detectivity (D*)	2.22 x 10 ¹⁴	2.00 x 10 ¹⁰	3.69 x 10 ¹⁴	6.67 x 10 ¹⁰
Bandwidth (-3dB – Hz, typical)	DC – 2 kHz	5 – 10 kHz	DC – 2 kHz	5 – 10 kHz
Power requirements	± 9 VDC to ± 15 VDC			
Connections	BNC signal output. Shielded power cable terminated with a DB-9 connector directly couples the unit with the PS/TC-1 Low Noise Power Supply / Controller.			


Mechanical Dimensions, Ambient and TE Housing

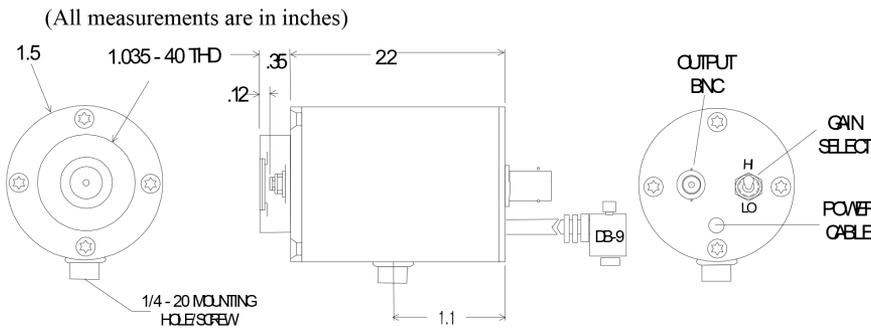

Electrical Diagrams, Ambient and TE Cooled

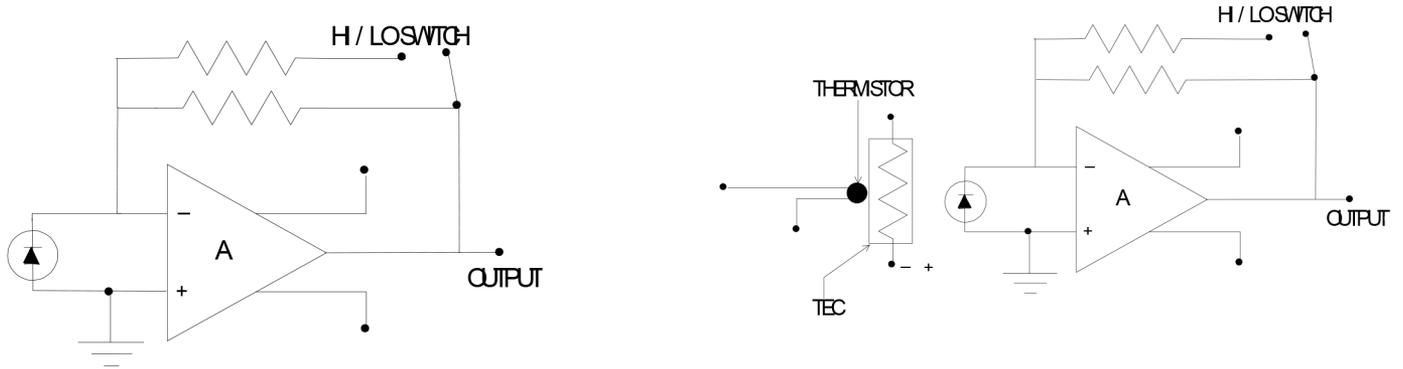

DB-9 Pin Out Diagrams, TE Cooled [Ambient]

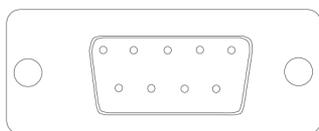

- | | |
|----------------------------|-------------|
| 1. Cooler (+) [No connect] | 6. +V |
| 2. Cooler (-) [No connect] | 7. -V |
| 3. Thermistor [No connect] | 8. GND |
| 4. Thermistor [No connect] | 9. Case GND |
| 5. No connect | |



OPTICAL BUILDING BLOCKS



HORIBA
Scientific

info.sci@horiba.com www.horiba.com/opticalbuildingblocks

USA: +1 732 494 8660
UK: +44 (0)20 8204 8142
China: +86 (0)21 6289 6060

France: +33 (0)1 69 74 72 00
Italy: +39 2 5760 3050
Brazil: +55 (0)11 5545 1500

Germany: +49 (0)89 4623 17-0
Japan: +81 (0)3 6206 4721
Other: +1 732 494 8660