


HORIBA

LabSpec 6

Spectroscopy Suite

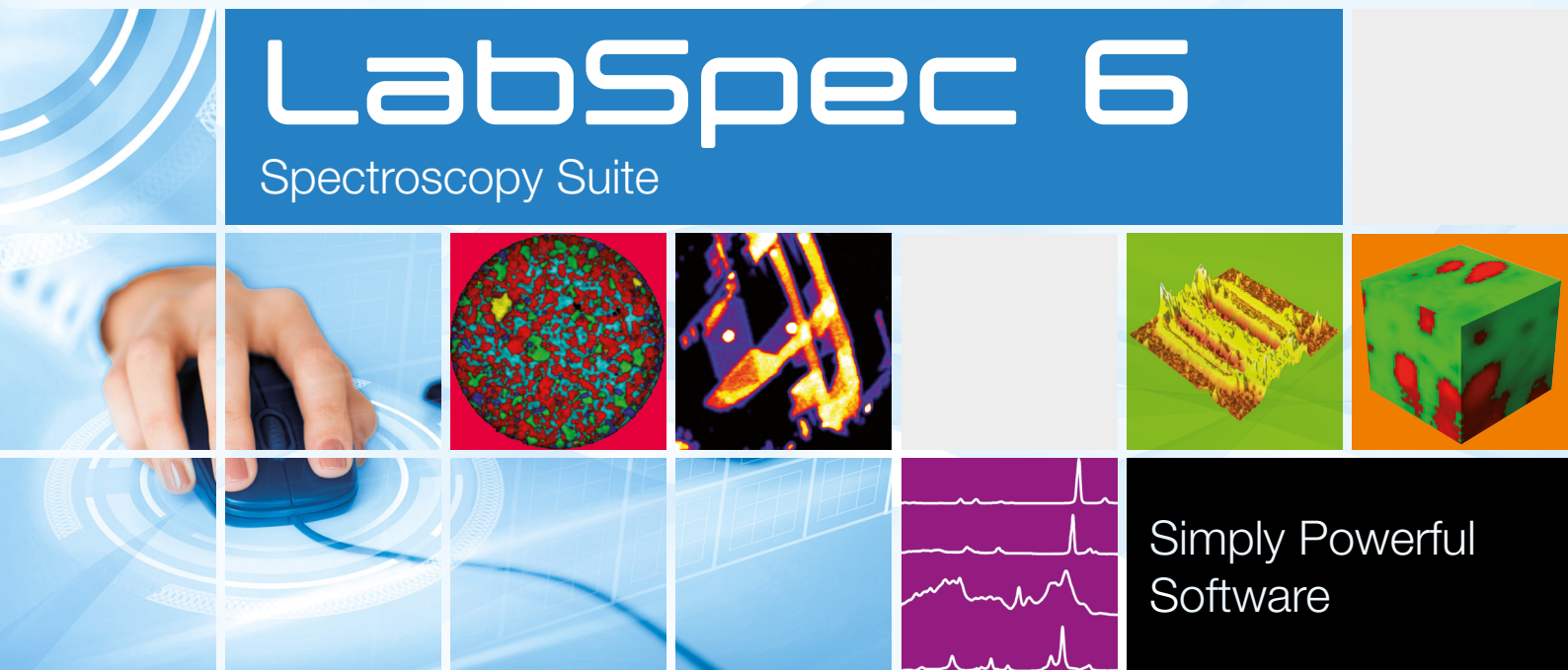

Simply Powerful
Software



LabSpec 6 Spectroscopy suite

The LabSpec 6 spectroscopy suite offers an intuitive platform for Raman, photoluminescence (PL) and cathodoluminescence (CL) spectroscopy and imaging.

The sixth generation of spectroscopy software from HORIBA Scientific builds on previous award winning platforms to present a unique experience in instrument control and data processing.

Within a single interface, LabSpec adapts to the needs of different users in multi-user lab environments, allows users to automate measurements / analysis and customize the user interface, and offers unrivalled functions to provide results fast.

LabSpec 6 software offers great modularity with exclusive LabStore Apps.

> Simplicity

An intuitive interface enables logical workflow through an experiment, from visualization of the sample and measurement set-up, to interpretation of the data, and final reporting of the results.

> Power

LabSpec is a powerful tool for the complete experiment, from instrument set-up and data acquisition, through to sophisticated routines for spectrum and image processing and display.

> Quality and Compliance

LabSpec 6 is validated software that contains all the tools for users to meet regulatory compliance including GxP and the FDA's 21 CFR part 11.


Instrument Control

LabSpec 6 is compatible with the full range of HORIBA Scientific's Raman spectrometers, including, analytical research LabRAM Soleil™, XploRA PLUS and the versatile research LabRAM HR Evolution microscopes. Their unique performance capabilities are made easily accessible via LabSpec's intuitive user interface.


- Full range, seamlessly glued **spectra** acquired from CCD, EMCCD, InGaAs and single channel (PMT, lock-in amplifier) detectors
- Automated **laser wavelength switching** with just a single mouse click ^{1,2,3}
- Auto-alignment: automatically check and **readjust the alignment** within a few seconds to compensate for changing environmental conditions ¹
- Optical **microscope imaging** with single field of view, and large area montage
- **SmartID™**: remove the risk of using the wrong objective or inappropriate acquisition parameters ¹
- Confocal **DuoScan™³ / QScan™¹** technology for adjustable laser spot size and high resolution sample mapping from deep UV to near infra-red
- Hyperspectral **surface and 3D volume mapping** using flexible point-by-point and ultra-fast SWIFT™ modes
- Automated multi-spectra experiment acquisition including **kinetic**, **polarization** and **heating/cooling** studies
- AFM-Server for combined **AFM-Raman** and **TERS** (Tip Enhanced Raman Spectroscopy) experiments
- **EasyNav™** Topographic and reflection **imaging** of the surface of tough samples

Workflows


LabSpec 6 software offers an intuitive interface which enables a logical workflow through an experiment, from visualization of the sample and measurement setup, to interpretation of the data, and final reporting of the results.

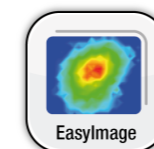

EasyImage ⁴ enables quick and easy acquisition of meaningful chemical images of your samples without needing a high level of Raman expertise

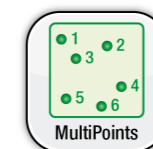

MultiPoints ⁴ provides automated Raman spectra acquisition. Create your own recipe by selecting multiple random positions on the sample

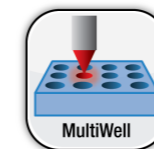

MultiWell ⁴ offers an easy interface for high throughput screening using traditional well plates, microtiter plates and other regularly arrayed sampling devices



ParticleFinder ⁴ offers a user-friendly tool for automated location, characterization and Raman analysis of particles. It is fully compatible with the latest market regulations

EasyNav™ Technology

With the groundbreaking EasyNav™ package, it is fast and easy to navigate in-focus, in real-time, to identify the region of interest and obtain sharp, clear Raman chemical images. HORIBA NavMap™ + NavSharp™ + ViewSharp™ apps can be used together or separately to deliver a powerful user experience for all Raman users.



SmartID automatically recognizes the microscope objective in use



NavMap helps to navigate around your large samples



NavSharp ⁴ provides real time focusing while navigating around samples




ViewSharp ⁴ constructs 3D sample topography and fully focused images



Mosaic enables fast imaging of large sample areas by video stitching

Spectral Acquisition

LabSpec 6 can acquire multi-dimensional datasets, full spectral range or custom ranges, including 1D profiles (such as line X, Y, depth Z, time, kinetic, temperature studies), 2D images (such as XY surface maps) and 3D cubes (i.e., confocal XYZ maps).

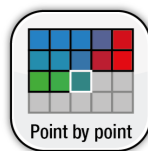

Point-by-point mapping provides ultimate sensitivity for imaging, ideal for low scattering materials



SWIFT mapping provides ultra-fast images with acquisition times below 500ms/point. Ideal for high scattering materials



SmartSampling ⁴ allows super large and super high resolution image acquisition up to 100 times faster, even for low scattering materials

Data Processing


Processing is a crucial step in spectroscopy. The preprocessing of raw spectral data helps to eliminate noise (and undesired signals,) along with the enhancement of discriminating spectral features. LabSpec 6 software includes a large palette of processing tools that can be applied in real time during acquisition or post-acquisition for enhanced qualitative and quantitative analysis!



FLAT enables Real-time on-the-fly fluorescence backgrounds removal from spectra



Data enhancement provides custom baseline, smoothing, normalization, thresholding and more

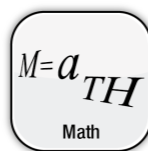

Math allows mathematical operations to be applied for spectral and image data manipulation


Analysis

Enhance the wealth of information by performing a complete Analysis of your spectral data!


LabSpec 6 guides you to Convert map data into chemical images!




chiSTain ⁴ is an innovative and intelligent application for fully automated processing and analysis of 2D Raman images

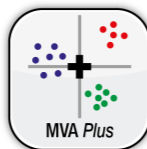

MVA Plus ⁴ combines unsupervised multivariate algorithms (PCA, MCR, HCA and K-means, CLS) for a quantitative and qualitative image generation

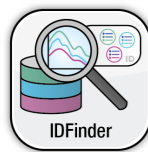

IDFinder provides spectral identification, mixture analysis, and spectral library management. It seamlessly integrates with all LabSpec6 applications, including ParticleFinder for precise particle identification

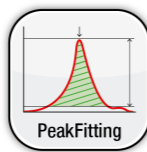

PeakFitting ⁴ analysis tool provides quantitative investigations of materials strain, interactions with the environment and crystallinity

Multimodalities


LabSpec 6 software is a flexible platform offering compatibility with multiple modalities.




AFM combines Raman and AFM systems, for push-button TERS and colocalized experiments



Epifluo provides compatibility with high performance, cooled fluorescence imaging cameras

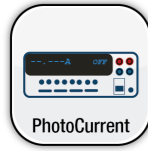

PhotoCurrent ⁴ allows the acquisition and display of photocurrent or photovoltage images

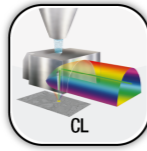

CL efficiently drives your SEM to achieve optimal colocalized cathodoluminescence analysis

App Packages


Save time and money by using a dedicated LabSpec Software application! An application package will combine all the processing and analysis tools that you need for a given application, thus automatically, in one click, you will get the answer to your question!



SI Stress ⁴ is a dedicated application for analysis of Silicon and photovoltaic cell samples



DLC ⁴ is a dedicated application for analysis of Diamond Like Carbon samples

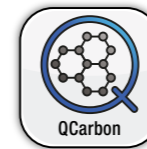

QCarbon simplifies Raman analysis for Carbon materials by processing spectral data and providing the ID/IG ratio.




Layers automates multi-layer material analysis, offering thickness and chemical identification for each layer.

Image Display

LabSpec 6 contains a wide variety of tools for working with hyperspectral Raman images and optical images. These include many advanced image processing tools and display functions for a perfect report!



DataOverlay combines all your multimodal sample images for hybrid sample analysis and viewing

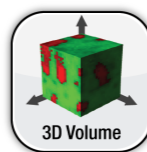

3D Volume integrated topographic surface and 3D confocal volume display with rotation and lighting controls



Image Enhancement offers a large palette of image processing tools for the enhancement of Raman and optical images

Reporting

Be ready to easily create optimized and customized measurement and analysis reports!



CustomReport customizes multi-data and multipage templates for fast reporting of results



Animation provides GIF animations of Raman and optical images for your presentation

Automation & Customization

Automate your own daily routines with LabSpec 6!



BatchProcess enables efficient fast batch data processing of multiple data files



OneClick allows fast, automated and easy Raman spectra acquisition



Custom GUI customizes your LabSpec 6 display to show just the tools you need in a multi-level user interface



Methods records routines allowing multiple measurements, processing and analysis steps to be fully and easily automated



Script & ActiveX allows in-app or external programming for instrument control and data processing

Compliance, Validation & Security

LabSpec 6 is a validated software designed to maximize ease of use, functionality and stability. It provides and ensure security and data integrity to enable compliant use of HORIBA Raman microscopes within regulated laboratory environments.



LabSpec 6 is a **validated software**, offering tested and documented functionality and performance



UserAccounts provides user access control and interface adaptation within multi-user environments



OneCheck ensure that your instrument is always perfectly aligned and calibrated



eVal validates Raman instruments for compliance with various pharmacopeias



ProtectionPlus ⁴ allows working in compliance with 21CFR Part 11 of the FDA's Code, with regard to electronic records and electronic signatures

Upgrades


Commercial LabSpec6 licenses include a free update plan to the latest version, through international service support. There are many benefits to keep your software up to date, including:

- Optimized user interface with clear, logical distribution of functions and controls; combined with the [User accounts](#) module it ensures even less-experienced users in large multi-user lab environments can utilize LabSpec's powerful functionality with ease.
- Access to the latest Apps such as [MVAPlus™](#) for multivariate analysis, [ParticleFinder™](#) for particle analysis, [χSTaiN™](#) for instant Raman analysis, [EasyNav™](#) for optimal sample navigation and much more.
- New and [improved capabilities](#) across the wide range of data acquisition, processing and analysis functions available in the LabSpec platforms.
- Create [stunning images](#) with a new selection of palettes and image display tools.
- [Faster data processing and display](#), utilizing LabSpec's multi-processing and multi-threaded core

Please contact your HORIBA Scientific service representative to enquire about upgrade requirements and costs for your system.

System Requirements	Recommended
CPU	Intel® Core™ i7
RAM	16 Gb or greater
Storage	SSD 100 Gb of free space
Monitor	1920 x 1080 or greater
Operating System	Windows® 10

Training and Support

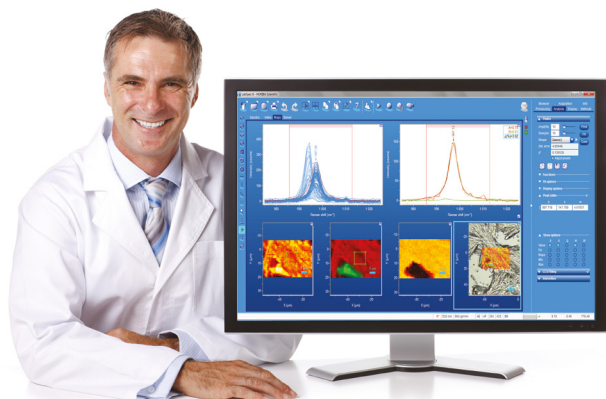

Our experienced Raman scientists based in well-equipped laboratories across the world are available to provide customized training both on-site and at HORIBA facilities.

Such training ensures confident and knowledgeable users who will be able to harness the full power of LabSpec 6 to obtain results with minimum effort.

- New user training
- Refresher courses
- Introduction to Raman
- "How-to" instruction
- Data interpretation
- Advanced techniques

Please contact your HORIBA Scientific representative for further information about available training.

Find out more at www.horiba.com/labspec6

HORIBA

info.sci@horiba.com

USA: +1 732 494 8660
UK: +44 (0)1604 542 500
China: +86 (0)21 6289 6060
Taiwan: +886 3 5600606

France: +33 (0)1 69 74 72 00
Italy: +39 06 51 59 22 1
India: +91 80 41273637
Brazil: +55 (0)11 2923 5400

www.horiba.com/scientific

Germany: +49 (0) 6251 8475 0
Japan: +81(75)313-8121
Singapore: +65 (0)6 745 8300
Other: +33 (0)1 69 74 72 00