HORIBA Global Application/Service Network

HORIBA





The HORIBA Group adopts IMS (Integrated Management System) which integrates Quality Management System ISO9001, Environmental Management System ISO14001, and Occupational Health and Safety Management System ISO45001. We have now integrated Business Continuity Management System ISO22301 in order to provide our products and services in a stable manner, even in emergencies.

Please read the operation manual before using this product to assure safe and proper handling of the product.

•The specifications, appearance or other aspects of products in this catalog are subject to change without notice

Please contact us with enquiries concerning further details on the products in this catalog.
Please contact us with enquiries concerning further details on the products in this catalog.
The color of the actual products may differ from the color pictured in this catalog due to printing limitations
It is strictly forbidden to copy the content of this catalog in part or in full.

The screen displays shown on products in this catalog have been inserted into the photographs through compositing

•All brand names, product names and service names in this catalog are trademarks or registered trademarks of their respective companies.

https://www.horiba.com/int/

HORIBA, Ltd.

Head Office 2 Miyanohigashi-cho, Kisshoin, Minami-ku, Kyoto.601-8510, Japan Phone: 81 (75) 313-8121 Fax: 81 (75) 321-5725

Tokvo Sales Office 2-6, KandaAwaji-cho, Chiyoda-ku, Tokyo, 101-0063, Japan Phone: 81 (3) 6206-4721 Fax: 81 (3) 6206-4730

HORIBA Taiwan, Inc. Taiwan

8F.-8, No.38, Taiyuan St. Zhubei City, Hsinchu County 30265, Taiwan (R.O.C.) Phone: 886 (3) 560-0606 Fax: 886 (3) 560-0550

HORIBA INSTRUMENTS (SHANGHAI) CO., LTD

China No.99, Chunxiu Rd, Anting Town, Jiading District, Shanghai, China 201804 Phone: 86 (21) 6952-2835 FAX: 86 (21) 6952-2823

HORIBA (China) Trading Co., Ltd. China Unit D. 1F. Building A. Synnex International Park, 1068 West Tianshan Road, 200335, Shanghai, China Phone: 86 (21) 6289-6060 Fax: 86 (21) 6289-5553 Beijing Branch 12F, Metropolis Tower, No.2, Haidian Dong 3 Street, Beijing, 100080, China Phone: 86 (10) 8567-9966 Fax: 86 (10) 8567-9066

HORIBA Instruments (Singapore) Pte Ltd. Singapore 3 Changi Business Park Vista #01-01, Singapore 486051 Phone: 65 (6) 745-8300 Fax: 65 (6) 745-8155

HORIBA Vietnam Company Limited Vietnam Lot 3 and 4, 16 Floor, Detech Tower II, No.107 Nguyen Phong Sac Street, Dich Vong Hau Ward, Cau Giay District, Hanoi, Vietnam Phone: 84 (24) 3795-8552 Fax: 84 (24) 3795-8553

PT HORIBA Indonesia Indonesia Japan Jl. Jalur Sutera Blok 20A, No.16-17, Kel. Kunciran, Kec. Pinang Tangerang-15144, Indonesia Phone: 62 (21) 3044-8525 Fax: 62 (21) 3044-8521 HORIBA KOREA Ltd.

> 25, 94-Gil, Iljik-Ro, Manan-Gu, Anyang-Si, Gyeonggi-Do, 13901, Korea Phone: 82 (31) 296-7911 Fax: 82 (31) 296-7913

HORIBA India Private Limited India 246, Okhla Industrial Estate, Phase 3 New Delhi-110020, India Phone: 91 (11) 4646-5000 Fax: 91 (11) 4646-5020 Technical Center D-255, Chakan MIDC Phase-II, Bhamboli Village, Pune-410501, India Phone: 91 (21) 3567-6000

HORIBA Instruments Incorporated 9755 Research Drive, Irvine, CA 92618, U.S.A. Phone: 1 (949) 250-4811 Fax: 1 (949) 250-0924

Ann Arbor Office 5900 Hines Drive, Ann Arbor, MI 48108, U.S.A. Phone: 1 (734) 213-6555 Fax: 1 (734) 213-6525 Sunnyvale Office 430 Indio Way, Sunnyvale CA94085, U.S.A. Phone: 1 (408) 730-4772 Fax: 1 (408) 730-8975 HORIBA New Jersey Optical Spectroscopy Center 20 Knightsbridge Rd, Piscataway, NJ 08854, U.S.A. Phone: 1 (732) 494-8660 Fax: 1 (732) 549-5125

HORIBA UK Limited

Kvoto Close Moulton Park Northampton NN3 6FL UK Phone: 44 (0) 1604 542500 Fax: 44 (0) 1604 542699

HORIBA (Austria) GmbH

Kaplanstrasse 5. A-3430 Tulln. Austria Phone: 43 (2272) 65225 Fax: 43 (2272) 65225-45

HORIBA FRANCE SAS

Korea

USA

UK

Austria

France 14 Boulevard Thomas Gobert - Passage Jobin Yvon CS 45002-91120 Palaiseau - France Phone: 33 (1) 69-74-72-00 Fax: 33 (1) 69-31-32-20

HORIBA ABX SAS France Parc Euromédecine, rue du Caducée, BP7290, 34184 Montpellier Cedex 4, France Phone: 33 (0) 4-67-14-15-16 Fax: 33 (0) 4-67-14-15-17

Germany

Sweden

talv

HORIBA Europe GmbH

Hans-Mess-Str.6, D-61440 Oberursel, Germany Phone: 49 (6172) 1396-0 Fax: 49 (6172) 1373-85 Leichlingen Office Julius-kronenberg-Str.9, D-42799 Leichlingen, Germany Phone: 49 (2175) 8978-0 Fax: 49 (2175) 8978-50

HORIBA Europe GmbH

Sweden Branch (Sodertalje) Sydhamnsvagen 55-57 SE-15138 Sodertalie Sweden Phone: 46 (8) 550-80701 Fax: 46 (8) 550-80567

HORIBA ITALIA SRI Torino Office Via Feroggio, 30, 10151 Torino, Italy

Phone: 39 (11) 904-0601 Fax: 39 (11) 900-0448 HORIBA Czech

Czech Prague Office Prumyslova 1306 / 7, CZ-10200, Praha 10, Czech Republic

Phone: 420 (2) 460-392-65 HORIBA 000 Russia

Altufievskoe shosse, 13, building 5, 127106, Moscow, Russia Phone: 7 (495) 221-87-71 Fax: 7 (495) 221-87-68

Laser Scattering Particle Size Distribution Analyzer



Bulletin:HRE-3683F

Laser Scattering Particle Size Distribution Analyzer





DCICIC LA-960V2 Series

Laser Scattering Particle Size Distribution Analyzer

HORIBA proudly introduces the newest breakthrough in particle size technology.

Laser Scattering Particle Size Distribution Analyzer Partica LA-960V2

This latest evolution in the LA series advances scientific knowledge for tomorrow's world through intuitive software, unique accessories, and high performance. The Partica LA-960V2 continues HORIBA's long standing tradition of leading the industry with innovative design in both the hardware and the software. The new optical design allows the user to visualize the particle dispersion in real time.



Proven high accuracy and resolution for wide application



Battery Capacitor 3D printing Paper Coating Pharmaceutical Cosmetics Food / Drink **Building materials**



Electrolyte

HORIBA's Original Optical Design

The groundbreaking optical design perfects the static light scattering particle sizing technique.

Advanced Detector Design

The number of detectors, angular range, and layout contribute to overall system performance.

The Partica LA-960V2 uses logarithmically spaced silicon photodiodes that detect a wide range up to 165 degrees to allowing the measurement of complete particle size distributions.

Automatic Laser Alignment in Seconds

Always make perfect measurements with computer-controlled laser alignment. The alignment process is completed in only a few seconds with HORIBA's innovative approach.









Superior Instrument-to-Instrument Precision

The Partica LA-960V2 is designed and built to provide the same experience regardless of manufacture date, operator skill, or geographic location. It achieves unmatched instrument agreement without the need for additional correlation procedures.

Guaranteed Accuracy

The Partica LA-960V2 is a capable of accurately measuring NIST-traceable size standards within 0.6% of specification. Fully compliant with ISO 13320 recommendations.

> Superior nanoparticle detection using red and blue light sources



0.05 µm 405 nm 0.07 µm 405 nm



0.05 µm 650 nm 0.07 µm 650 nm

DCITICCI LA-960V2 Series

Laser Scattering Particle Size Distribution Analyzer

State of the Art Sampling Systems

Wet Measurement

Sample-to-sample analysis in less than 60 seconds

The Partica LA-960V2 wet circulation system is an easy, fast and very powerful dispersion system. The standard wet system offers a full package which includes dispersant fill pump, liquid level sensor, circulation pump, 30 W in-line ultrasonic probe, and drain valve. It is all software-controlled for true one-button operation. This advanced design provides highly reproducible particle size results.





Fluid sensor

Dry Measurement

Automated, powerful dry powder dispersion

The Partica LA-960V2 Powderjet combines several unique and patented features to provide the most reproducible dry measurements. Use the Auto Measurement function to control vacuum, air pressure, powder flow, start/stop conditions, measurement duration, and data processing. Designed to handle every application including small sample amounts, friable powders, and highly agglomerated materials.

Smart Scans - Trigger Function

This function allows precious materials to be measured accurately. When the sample amount is limited or has low flowability, the trigger functions starts and stops the measurement with faultless precision.

Self-adjusting Powder Flow

Historically, the biggest challenge in dry powder measurement involved establishing an even powder flow. The Partica LA-960V2 Powderjet has solved that challenge with a self-adjusting feedback loop to maintain a constant laser transmittance. This is a crucial factor in creating reliable, reproducible dry powder size measurements

Available chutes

Standard chute Equipped with every Powderjet Dry Feeder and used for most powders.



V type chute

This chute is ideal for both small amounts of powder and powder which does not flow easily, e.g. magnetic powders.

Compressed air Air guidance plates Cell window Laser light Vacuum

Sample

Coated chute

This coated chute is useful for samples which adhere to the stainless steel standard chute

Accessories

High Concentration Cell

Feature



The high concentration cell unit allows measurement closer to the original concentration with low dilution rates, no dilution and variable concentration.

Typical Applications

Understand the particle dispersion state of high concentration slurry, such as positive and negative electrode materials of batteries, inks. paints, pigments, emulsions etc.

Feature

Imaging Analysis Unit (Built-in option)



The imaging unit visualizes the particles in the wet circulation system. It is very small and built into the main unit, and can acquire particle images information without changing the usability.

Measurement range : 9 µm-1000 µm

MiniFlow (Circulation system)

Feature



The MiniFlow minimizes sample and dispersant amounts This miniaturized circulation system features fill and circulation pumps, an ultrasonic probe, and drain valve for fully automated operation. Measurement range: 10 nm - 1000 µm

Typical Applications

Valuable samples requiring powerful dispersion and materials requiring hazardous dispersants such as organic solvent and oil solvent

Auto Sampler (Wet measurement in powder)



The Auto Sampler is a rotary table-type automatic sampling system equipped with 24 detachable sample cups.

Vacuum sampler This accessory is useful for measuring very small amounts of powder. *Sampling table is included



Comparison of the results of undiluted suspension measurement and flow measurement (diluted with dispersant)



Typical Applications

Find and count small amounts of large particles and aggregated particles. 8 types of particle shape analysis and the alarm of the observed "bubble" are also possible

Fraction Cell



Feature

The Fraction Cell makes measurements with only micrograms of sample. This unique accessory is available in 5, 10 and 15 mL volumes and fully solvent resistant

Typical Applications

Samples requiring minimal dispersion such as precious samples, bio material and highly volatile solvent. Samples to measure without dispersing force.



Laser Scattering Particle Size Distribution Analyzer

Innovations in hardware and software

Performance

State of the art nanoparticle measurement

The advanced design of the Partica LA-960V2 allows for easy measurement of nanoparticle applications. NIST-traceable size standards verify that the Partica LA-960V2 accurately measures peaks as fine as 30 nanometers.



Overlay of 30, 40, 50 and 70 nm results

Speed One click measurement Navigation system

This incredible speed is made possible by automatic laser alignment, fully automated liquid handling and intuitive software design.



Wide range

Measurement range 10 nm - 5000 µm

The Partica LA-960V2 features a wide measurement range to measure every application. The unique optical bench is user-friendly and standard in every Partica LA-960V2 configuration.

Operation Intuitive software

The Partica LA-960V2 software is designed to be intuitive, allowing the user to check particle size distribution in real time. There is also a navigation system that allows one click measurements.

In addition, Method Expert software makes it easy to create robust, powerful methods for research and development purposes as well as quality control. The Method Expert is a series of guided, automated tests with advice to help the user choose values for refractive index, concentration, ultrasonic dispersion, pump speed and measurement duration. Without any training, users can generate effective data in a short amount of time using the software.



Measurement Screen Real-time display of the measurement result



The Method Expert recommends the most suitable refactive index

Data support

Traceability certification

- 21 CFR Part 11 compliant support optional software
- ✓ IQ/OQ/PQ documents support
- Data correlation support with old model

Partica LA-960V2 Standard Model

Measurement Principle	Mie scattering and Fraunhofer diffraction
Measurement Range	10 nm - 3000 μm
Measurement Time	Typical measurement takes 60 seconds from liquid filling, sampling and measurement to rinsing
Measurement Method	Circulation measurement or fraction cell measurement (Fraction cell is optional)
Sample Quantity	Approximately 10 mg - 5 g (Depending on the particle size, distribution and density)
Dispersing Volume	Approximately 180 mL for standard pumping system 5/10/15 mL for Fraction Cell accessory Minimum volume 35 mL for Mini Flow accessory Approximately 1 L of LiterFlow option
Available Carrier Fluid	Aqua* (A type), Organic solvent (S type) (*Small volume ethanol can be used as a dispersing additive)
Communication	USB 2.0
Light Sources	Red solid state 5 mW laser diode (650 nm), Blue solid state 3 mW LED (405 nm)
Dispersion System	In-line ultrasonic probe: 30 W, 20 kHz, adjustable levels Circulation pump: Fully automated fill and circulation pumps, 4 selectable fill levels, 15 selectable circulation speeds (max: 10 L/min)
Operating Conditions	15 - 35°C (59 - 95°F), relative humidity 85% or less (No condensation)
Power	AC100-240 V 50/60 Hz, 300 VA
Dimensions	705 (W) x 565 (D) x 500 (H) mm
Mass	54 kg
Computer Requirements	PC operation, Software compatible with Windows® 10 and 11 64-bit environments, *Contact HORIBA for additional operating system compatibilities

With Powderjet Dry Feeder Accessory

orating feeder cuum-driven evacuation
cuum-driven evacuation
0 nm - 5000 μm
ommunication: Serial cable to the main unit easurement: Vibrating feeder controlled automatically via feedback or manually by user, Vacuum AUTO/OFF, Compressed Air AUTO/OFF, Air pressure adjustable from 0 - 0.4 MPa in 40 steps
pical measurement takes 2 seconds or longer
- 35°C (59 - 95°F), ative humidity 85% or less (no condensation)
2 (W) x 321 (D) x 244 (H) mm ot including dimensions of projections and the measurement unit)
C100 V, 120 V, 230 V, 50 or 60 Hz, 1500 VA cluding vacuum but the measurement unit)
mpressed air supply origin pressure: 0.4 - 0.8 MPa mpressed air controlling range: 0.01 - 0.4 MPa
ick connector for resin tube with 6 mm outer diameter ompressed air supply equipment must be provided separately)
cuum is equipped as standard

* Manufacturers and models indicated for vacuum, air compressor, computer, monitor and/or printer are subject to change

Air Compressor Inlet pressure within 0.5 - 0.8 MPa, Tank capacity 26 L or larger, Flow rate 45 L/min or faster

Dimensions (mm)



Class 1 Laser Product

Dimensions (mm)



Class 1 Laser Product