

www.horiba.com/us/particle

Jeffrey Bodycomb, Ph.D.

Introduction to Dynamic Light Scattering for Particle Size Determination



© 2016 HORIBA, Ltd. All rights reserved Explore the future

Automotive Test Systems | Process & Environmental | Medical | Semiconductor | Scientific

HORIBA



Sizing Techniques





Two Approaches to Image Analysis

Dynamic: particles flow past camera



1 – 3000 um

Static: particles fixed on slide, stage moves slide



0.5 – 1000 um 2000 um w/1.25 objective

© 2016 HORIBA, Ltd. All rights reserved.



Laser Diffraction



Laser Diffraction

- •Particle size $0.01 3000 \ \mu m$
- •Converts scattered light to particle size distribution
- Quick, repeatable
- Most common technique
- •Suspensions & powders

© 2016 HORIBA, Ltd. All rights reserved.



HORIBA



Laser Diffraction



HORIBA



What is Dynamic Light Scattering?

- Dynamic light scattering refers to measurement and interpretation of light scattering data on a <u>microsecond</u> time scale.
- Dynamic light scattering can be used to determine
 - Particle/molecular size
 - Size distribution
 - Relaxations in complex fluids



Other Light Scattering Techniques

- Static Light Scattering: over a duration of ~1 second. Used for determining particle size (diameters greater than 10 nm), polymer molecular weight, 2nd virial coefficient, R_g.
- Electrophoretic Light Scattering: use Doppler shift in scattered light to probe motion of particles due to an applied electric field. Used for determining electrophoretic mobility, zeta potential.



DLS Optics



HORIBA



Brownian Motion

Particles in suspension undergo Brownian motion (random thermal motion.





- Brownian Motion
 - Random
 - Related to Size
 - Related to viscosity
 - Related to temperature



DLS signal

 Random motion of particles leads to random fluctuations in signal (due to changing constructive/destructive interference of scattered light.



time (microseconds)



Correlation Function

 Random fluctuations are interpreted in terms of the autocorrelation
 2.0 γ
 function (ACF), C(τ).



 $C(\tau) = 1 + \beta \exp(-2\Gamma \tau)$



Gamma to Size



13 HORIBA



What is Hydrodynamic Size?

 DLS gives the diameter of a sphere that moves (diffuses) the same way as your sample.



HORIBA



Hydrodynamic Size

- The instrument reports the size of sphere that moves (diffuses) like your particle.
- This size will include any stabilizers bound to the molecule (even if they are not seen by TEM).

Gold Colloids

Technique	Size nm
Atomic Force Microscopy	8.5 ± 0.3
Scanning Electron Microscopy	9.9 ± 0.1
Transmission Electron Microscopy	8.9 ± 0.1
Dynamic Light Scattering	13.5 ± 0.1

SEM (above) and TEM (below) images for RM 8011







Nanogold Data







Nanogold Data



© 2016 HORIBA, Ltd. All rights reserved.

Explore the future



Lab to Lab comparison

Colloidal Silica

	Mean determined Z-average size (nm)	COV (%)
Dynamic Light Scattering with SZ-100, laboratory 1	34.4	0.7
Dynamic Light Scattering with SZ-100, laboratory 2	34.6	0.3

HORIBA



Emulsion Polymers

IUPAC definition Emulsion polymerization:

Polymerization whereby monomer(s), initiator, dispersion medium, and possibly colloid stabilizer constitute initially an inhomogeneous system resulting in particles of colloidal dimensions containing the formed polymer.





Polystyrene Latex Sample

	Measurement <u>M</u> anual Measurement	Loois Heip		
Size •	<u> </u>			
Particle Size	Data	Comple Nome	Contraring Angle	Z Average(pm) DI
Legena No. D.Q.C	Date Thursday, May 26, 2011 12:12:10.1	Sample Name	Scattering Angle	2-Average(nm) PI
	Hitistiay, May 20, 2011-12, 12, 10	-MIPSE 100 mm 4 mmsodeg	90	105.4 0.012
Average				105.4 0.012
8. D.				
Circ Diskhafen I. e	la col			
Size Distribution Autocorrelation Functio	n Hesidual	Condition Table		
20	-100	Measurement Res	ults	^
20-		Sample Name	: PSL 100	nm 4 min90deg
		ID#	: 2011052	60912010
~ 15	N	Measurement Type	: Inurso: Particle	ay, May 26, 2011 12:12
op 15		ND Filter 1	: No.8 3%	T
~		Temperature of the Hold	er : 25.2 °C	
~	- .	Particle Name	: mono-p	olystyrene
Cod 10		Divis Division di a sa	· 240 Cod	onds
0 4 9 10	-40	Run Duration	. 240 Sec	an an I. I. and a data for the state of the
10	-40 [%] ap	Run Duration Representation of Result	: Scatteri	ng Light Intensity
Го иел 10- ль з ль з ль	-40 р - Ц - Ц	Representation of Result Calculation Result	: 240 Sec : Scatteri S	ng Light Intensity
Со и 10- в - и	-40 ^H -40 - H 20	Calculation Result Cumulant Operation	S	ng Light Intensity
Le dre Domen de Le composition de la composition	-40 ^H app - D - 20	Representation of Result Calculation Result Cumulant Operation Z-Average	: 240 Sec : Scatteri S : 105.4 ni	ng Light Intensity m
To the second se	-40 ⁹ 9 - ¹ 9 - 20 20	Representation of Result Calculation Result Cumulant Operation Z-Average Pl	: 240 Sec : Scatteri S : 105.4 ni : 0.012	ng Light Intensity m
10 10 10 10 10 10 10 10 10 10	-40 - 9 - 9 - 9 - 20 - 20 - 1000 10000	Representation of Result Calculation Result Cumulant Operation Z-Average PI Cumulants S. D.	: 240 Sec : Scatteri S : 105.4 ni : 0.012 : 11.5 nm	ng Light Intensity m
10 I I I I I I I I I I I I I I I I I I I	-40 	Representation of Result Calculation Result Cumulant Operation Z-Average PI Cumulants S. D.	: 240 Sec : Scatteri S : 105.4 ni : 0.012 : 11.5 nm	ng Light Intensity m
10 10 10 10 10 10 10 10 10 10 10 10 10 1	-40 99 - 99 - 20 20 100 10000	Representation of Result Calculation Result Cumulant Operation Z-Average PI Cumulants S. D.	: 240 Sec : Scatteri S : 105.4 ni : 0.012 : 11.5 nm	ng Light Intensity m

HORIBA



Polydisperse Sample Cumulants

For a mixture of sizes, the autocorrelation function can be interpreted in terms of cumulants. This is the most robust method of analyzing DLS data.

$$C(\tau) = 1 + \beta \exp(-2\Gamma\tau)$$

$$C(\tau) = 1 + \beta \exp\left[2\left(-\overline{\Gamma}\tau + \left(\frac{\mu_2}{2!}\right)\tau^2 - \cdots\right)\right]$$

$$\overline{\Gamma} = \overline{D_m}q^2 \qquad \text{"z-average size"}$$

$$D_{z,h} = \frac{k_B T}{3\pi\eta(T)\overline{D_m}} \qquad Polydispersity = \frac{\mu_2}{\Gamma^2}$$

$$e^{2t} = 2t^2 + 2t^2 + 2t^2 + 2t^2 + 2t^2$$



Run	Z-average Diameter (nm)	Polydispersity Index
1	473.2	0.127
2	479.5	0.066
3	478.8	0.077
4	487.7	0.039
Avg.	479.8	0.077

22 HORIBA



Mixtures of Particles



Sum the autocorrelation functions

23 HORIBA



Polydisperse Sample (ILT)

A more general relationship can be given between the autocorrelation function and the size distribution. Let each size have a relation constant Γ . The scattering from each population is then given by S(Γ). Now we have an integral equation. Solving for S(Γ) gives us size distribution.

$$C(\tau) = 1 + \beta |g^{(1)}(\tau)|^2$$





Bimodal Sample

Nominal 20 nm and 500 nm latex run individually





Bimodal Sample

Mixed sample (in black)







Dust

- Dust: large, rare particles in the sample
- Generally not really part of the sample
- Since they are rare cannot get good statistics



Filtering

- Filter to remove dust. If particles are too large (D >50 nm for 0.1 μm filter), at least filter diluent.
- Filters available in sizes 20nm to 2μm
- We can also centrifuge the sample and extract the supernatant.



© 2016 HORIBA, Ltd. All rights reserved.



Settling and DLS

Particle Diameter (μm)	Movement due to Brownian Motion		Movement due to Gravitational Settling
0.01	2.36	>>	0.005
0.25	1.49	>	0.0346
0.50	1.052	>	0.1384
1.0	0.745	~	0.554
2.5	0.334	<	13.84
10.0	0.236	<<	55.4

The Natural limit for Dynamic Light Scattering: Gravitational Settling

Gravitational Settling occurs at about 1-3µm

Explore the future



Why DLS?

- <u>Non-invasive</u> measurement
- Requires only <u>small quantities</u> of sample
- Good for <u>detecting trace amounts</u> of aggregate
- Good technique for <u>macro-</u> <u>molecular sizing</u>



Nanoparticle Analyzer

 Single compact unit that performs size, zeta potential, and molecular weight measurements.



© 2016 HORIBA, Ltd. All rights reserved.

Explore the future

Automotive Test Systems | Process & Environmental | Medical | Semiconductor | Scientific

31 HORIBA



Ask a question at <u>labinfo@horiba.com</u>

Keep reading the monthly <u>HORIBA Particle</u> e-mail newsletter!

Visit the **Download Center** to find the video and slides from this webinar.

Jeff Bodycomb, Ph.D.

- P: 800-446-7422
- E: jeff.bodycomb@horiba.com

