

Password Reset Procedure for LAQUA 1000 Series Bench Meters

If you've forgotten the password for your PH1100, PH1200, PH1300, EC1100, or PC1100 meter, follow these steps to reset the password.

1 Press the SET key and MODE key simultaneously for 5 seconds.




2 Press ENT key.



3 Press SET key.



4 Press down arrow key until GEN appears on the screen.

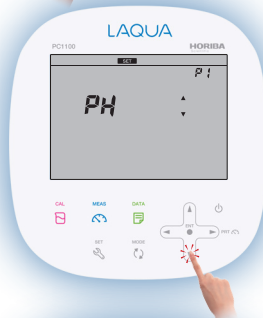

5 Press ENT key.




6 Press down arrow key again until PASS appears on the screen.



7 Press ENT key twice.



8 Enter a new 4-digit password using the arrow keys: left/right arrows—to move to the left or right digit, up/down arrows—to increase or decrease the digit.

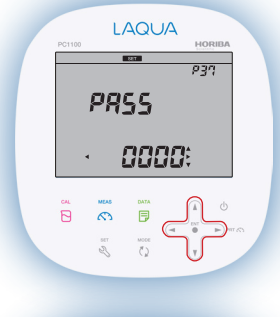

9 Press ENT key to save the new 4-digit password.



10 New password is set.

